

## SPECIFICATIONS

### Range:

–58 to 302°F (–50 to 150°C)

### Resolution

0.1° from –20.0 to 100°C

### Accuracy:

±1°C from –20 to 200°C

## OPERATION

1. Slide the power “ON/OFF” switch to the “ON” position.
2. Slide the “C/F” switch to select the desired reading.
3. Press “HOLD” button once to freeze the display.
4. Press the “HOLD” button again to resume measurement.
5. Slide the power “ON/OFF” switch to the “OFF” position to conserve battery life.
6. Keep plastic housing and plastic handle away from heat source. Do not place inside oven or expose to open flame.
7. Do not immerse probe handle or cable in liquid.

## BATTERY REPLACEMENT

1. Use a coin to open the battery cover on back of unit.
2. Replace with catalog number 1039.

**WARRANTY, SERVICE, OR  
RECALIBRATION**

For warranty, service, or recalibration, contact:

**CONTROL COMPANY**

12554 Old Galveston Rd. Suite B230

Webster, Texas 77598 USA


Ph. 281 482-1714 • Fax 281 482-9448

E-mail [sales@control3.com](mailto:sales@control3.com)

[www.traceable.com](http://www.traceable.com)

Control Company is ISO 9001:2008 Quality-Certified  
by DNV and ISO/IEC 17025:2005 accredited as a  
Calibration Laboratory by A2LA.

Traceable® is a registered trademark of  Control Company.

©2016  Control Company. 92-4146-00 Rev. 1 061416

**TRACEABLE®  
MINI-  
THERMOMETER  
INSTRUCTIONS**